

# WF-10 Dino-Lite Connect

Dino-Lite Connect is a wireless device designed for Dino-Lite connected to your smartphone or tablet with iOS or Android OS.



## Summary

Dino-Lite Connect is released as an innovative alternative to conduct multiple tasks besides working with wired products. It is an option allowing smartphones or tablets to transmit or receive data using WiFi.

Specializing in being utilized with smartphones or tablets, its app, DinoConnect featuring multi-functions, provides you a free-and-easy way to supervise real-time status instead of staying on a designated workstation. In addition, the remote control gives you the capability executing any command for Dino-Lite that is away from you.

### **Note :**


System Requirement  
Android 4.1.1 or later  
iOS 6.x or later

The Dino-Lite models compatible to the WF-10 and DinoConnect are the Dino-Lite Premier (R4) series and Edge series and with USB interface

## Gallery

[Sample Photos]


## Package Contents

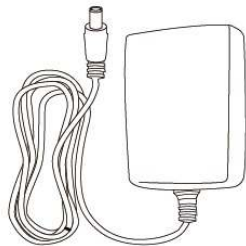

**Adaptor**




**WF-10 Dino-Lite Connect**

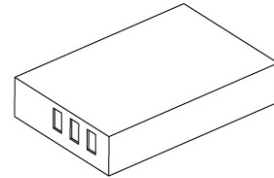




**Battery**

## Overview

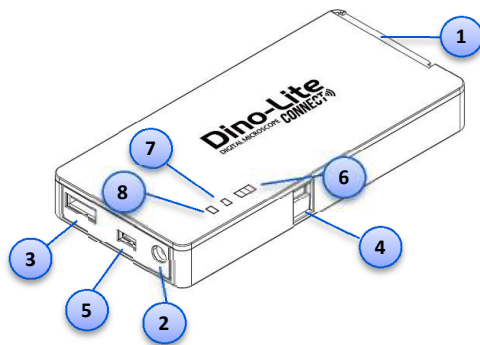

- |                           |                               |
|---------------------------|-------------------------------|
| 1. Battery Slot           | 5. Mini USB Port for Charging |
| 2. DC Input for Charging  | 6. Power Indicator            |
| 3. USB Port for Dino-Lite | 7. WiFi Indicator             |
| 4. Switch On/Off          | 8. Switch Indicator           |

## Hardware Preparation

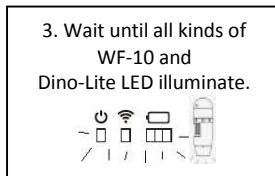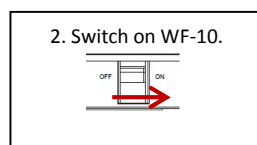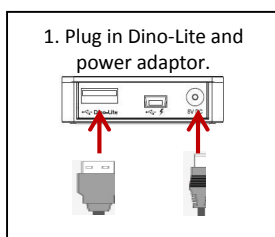


**Note :** The mini USB cable is excluded from our package contents. The Dino-Lite models compatible to the WF-10 and DinoConnect are the Edge series. Using supplied battery and adaptor is recommended.

## Using Dino-Lite Connect

**For Portable usage :**

After performing the said steps in Hardware Preparation, also perform the following

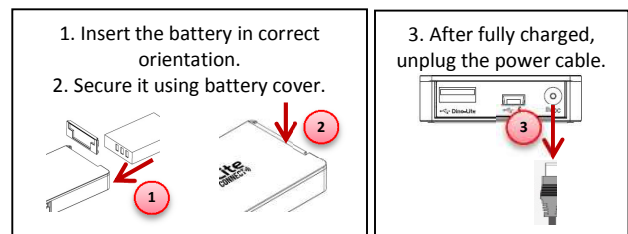

**Note :** Make sure that the battery is charged for at least 8 hours prior to initial use.

**For stationary usage :** Battery is not a strictly necessary item. You may perform the said step 1. through 3. in Hardware Preparation only, whereas it is still functional.

## DinoConnect Application Download

- To download DinoConnect, visit App Store for iOS version or Google Play for Android version.
- Search for free "DinoConnect" app and download.
- Select "Settings" on your wireless devices, then select "Wi-Fi"
- A list of available networks appears. Select the default "Dino-Lite WiFi Streamer" (The SSID is editable using DinoConnect App)
- Run DinoConnect App.