

## Silent Installation and Uninstallation Instructions

---

Instructions on how to create a silent installation and uninstallation file for DinoCapture 2.0 software or device driver of the Dino-Lite Pro series.

Notice: The AM211, AM311, and AM313 series may not be suitable for silent installation.

### Silent Installation

#### Recording the response File

1. Download the DNC\_SI from: [http://www.dino-lite.com/support/DNC2\\_SI\\_v1528A.exe](http://www.dino-lite.com/support/DNC2_SI_v1528A.exe)
2. Extract the DNC2\_SI.exe to the desired directory (C:\ is recommended).
3. Access the command prompt (cmd) from Windows.
4. Change to the directory where contains the setup scrip.

For example: cd C:\DNC2\_SI\_v1525A\DNC2\_AP

```
C:\Users\ANMO>cd C:\DNC2_SI_v1525A\DNC2_AP
```

5. Run the installer with switch /r (setup.exe /r).  

```
C:\DNC2_SI_v1525A\DNC2_AP>setup.exe /r
```
6. Proceed through the dialogs and complete the installation. This action creates a setup.iss response file under C:\Windows.
7. Put the setup.iss into the DNC2\_AP file.

#### Play back a response file

Copy the DNC2\_AP (with setup.iss) file to the target computer and run the installer through command prompt with the /s command (setup.exe /s). This will perform an unattended installation.

```
C:\DNC2_SI_v1525A\DNC2_AP>setup.exe /s
```

Note: If plan to create a silent response file for the Dino-Lite Pro series device driver, repeat steps 3 to 6 but change the directory to DN4xxDrv.

### Silent Uninstallation

#### Recording the response file

1. Access the command prompt (cmd) from Windows.
2. Change to the directory which contains the setup scrip.

For example: cd C:\DNC2\_SI\_v1525A\DNC2\_AP

```
C:\Users\ANMO>cd C:\DNC2_SI_v1525A\DNC2_AP
```

3. Run the uninstaller with setup /r /f1 (with the directory path where to create

uninstall.iss). Proceed through the dialogs and complete the uninstallation. This action creates a “uninstall.iss” response file in the designated location.

For example: setup /r /f1C:\DNC2\_SI\_v1525A\DNC2\_AP\uninstall.iss

```
C:\DNC2_SI_v1525A\DNC2_AP>setup /r /f1C:\DNC2_SI_v1525A\DNC2_AP\uninstall.iss
```

#### Play back a response file

Copy the DNC2\_AP (with uninstall.iss) to the target computer, and run the uninstaller through command prompt. This will perform a silent uninstallation.

For example: setup /s /f1C:\DNC2\_SI\_V1525A\DNC2\_AP\uninstall.iss

```
C:\DNC2_SI_v1525A\DNC2_AP>setup /s /f1C:\DNC2_SI_U1525A\DNC2_AP\uninstall.iss
```

Note: If plan to create a silent response file for the Dino-Lite Pro series device driver, repeat steps 1 to 3 but change the directory to DN4xxDrv.