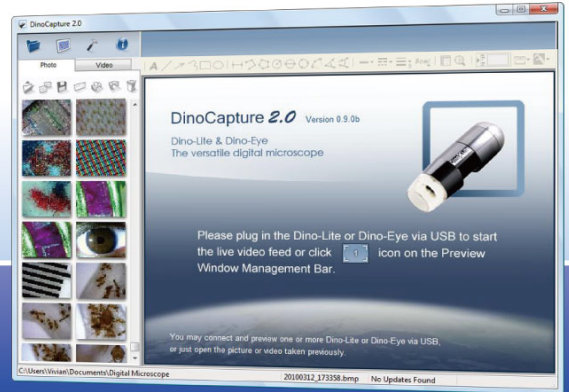
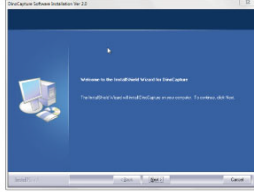

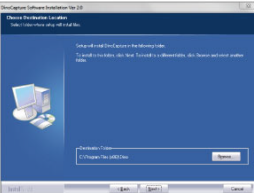
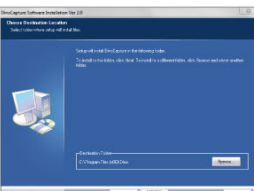
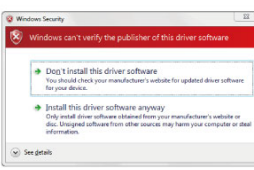
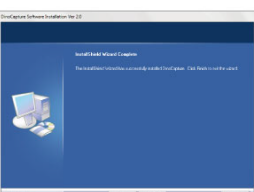


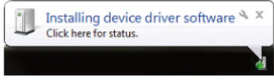
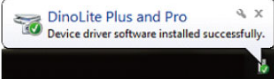
# DinoCapture 2.0

## Quick Installation Guide



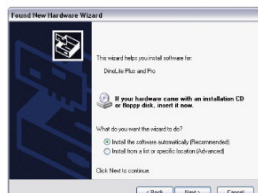
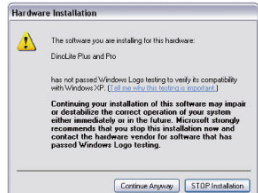
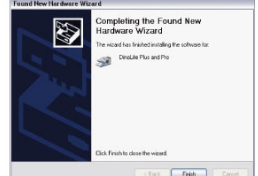
<http://www.dino-lite.com>



- 1 Power on your PC and insert DinoCapture 2.0 installation CD into CD-ROM drive. The installation wizard should start automatically.
 
- 2 Click "Next" and the Install Shield Wizard for the DinoCapture 2.0 will start. (The "Open File-Security warning" may appear on some systems. Select "Run")
 
- 3 Read the User License Agreement. If agree, press "Yes" to continue or "No" to stop the installation.
 
- 4 Select a destination folder for the DinoCapture 2.0 software. When done press "Next"
 
- 5 Click "Install" to start installing the software.
 
- 6 If the Windows Security warning message appears, click "Install this driver software anyway".
 
- 7 When the installation is complete, selecting "Finish" completes the software installation.
 

- 8 Now plug in the Dino-Lite to start the Driver Installation. A bubble will pop up saying "Installing device driver software". **Please WAIT** until Windows says "Device driver software installed successfully" for completed device driver installation.
 

- 9 Start DinoCapture 2.0 by double clicking the desktop icon.\*
 

**\*Windows XP Users continue reading steps 10-15 for Driver Software Installation instructions-Found New Hardware Wizard.**

- 10 Plug in the Dino-Lite to start the Found New Hardware Wizard.
 
- 11 Select the "No, not this time" radial button. And then press "Next".
 
- 12 Select "Install the software automatically" radial button then press "Next"
 
- 13 If the Windows Logo testing warning appears select "Continue Anyway".
 
- 14 Click "Finish" to complete the installation.
 
- 15 You can now use the Dino-Lite with the DinoCapture software.

Note: Please refer to the User's Guide in the DinoCapture 2.0 "Help" menu to learn more about DinoCapture 2.0 software.